## ***Installation Guide***

***Guida d'installazione / Guide d'installation***

***Guía de instalación / Installationsanleitung***

**UNIVERSAL H-STAND**

**30586771**

**TELESCOPIC H-STAND**

**30586772**

## General Information / Informazioni Generali / Informations Générales / Información General / Allgemeine Hinweise

Before using the unit, please read the following instruction manual carefully.  
Prima dell'utilizzo dello strumento si raccomanda di leggere attentamente il seguente manuale operativo.  
Avant d'utiliser l'instrument, il est recommandé de lire attentivement le présent manuel d'instructions.  
Antes de utilizar el instrumento, le recomendamos que lea con atención el siguiente manual de funcionamiento.  
Bitte lesen Sie vor Inbetriebnahme des Geräts diese Bedienungsanleitung sorgfältig durch

**This unit must be used for laboratory applications indoor only.** The manufacturer declines all responsibility for any use of the unit that does not comply with these instructions. If the product is used in a not specified way by the manufacturer or with not specified accessories, product's safety may be compromised.

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**This unit has been designed and manufactured in compliance with the following standards:**

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The adjustment of the screw can be used to compensate for unevenness on the set up plate; use the lock nut to lock the adjustment screw.

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Check the mounting position of the stirrer to avoid the danger of the unit tipping over.

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Observe the maximum permissible load-bearing capacity ( see "Technical features").

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Follow the safety instructions contained in the operating instructions for the stirrer.

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Before first use and periodically verify that the instrument is securely attached to the stand with the clamp and the beaker with the sample are secure. Vibration generated during general use may cause loosening.

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In relation to multiple applications and configurations, it's not possible to guarantee the user safety on the base of construction requirements established for the product. Therefore additional safety measures may be necessary for the user in relation to the specific application.

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Pay attention to the screw handle for stand height lock (Pos 3 Figure N°1)

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Pay attention during the height adjustment of the stand mobile part regulation making sure to not interfere with devices fixed to the support and other objects connected to the process for example limit switch vessel etc.

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Pay attention when the height lock handle is unscrewed: the stand mobile shaft has to be held and secured with the other hand to avoid it moving unexpectedly.

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Always store the stand in a relaxed state.

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The telescopic tube must not be directed at people, objects and/or animals.

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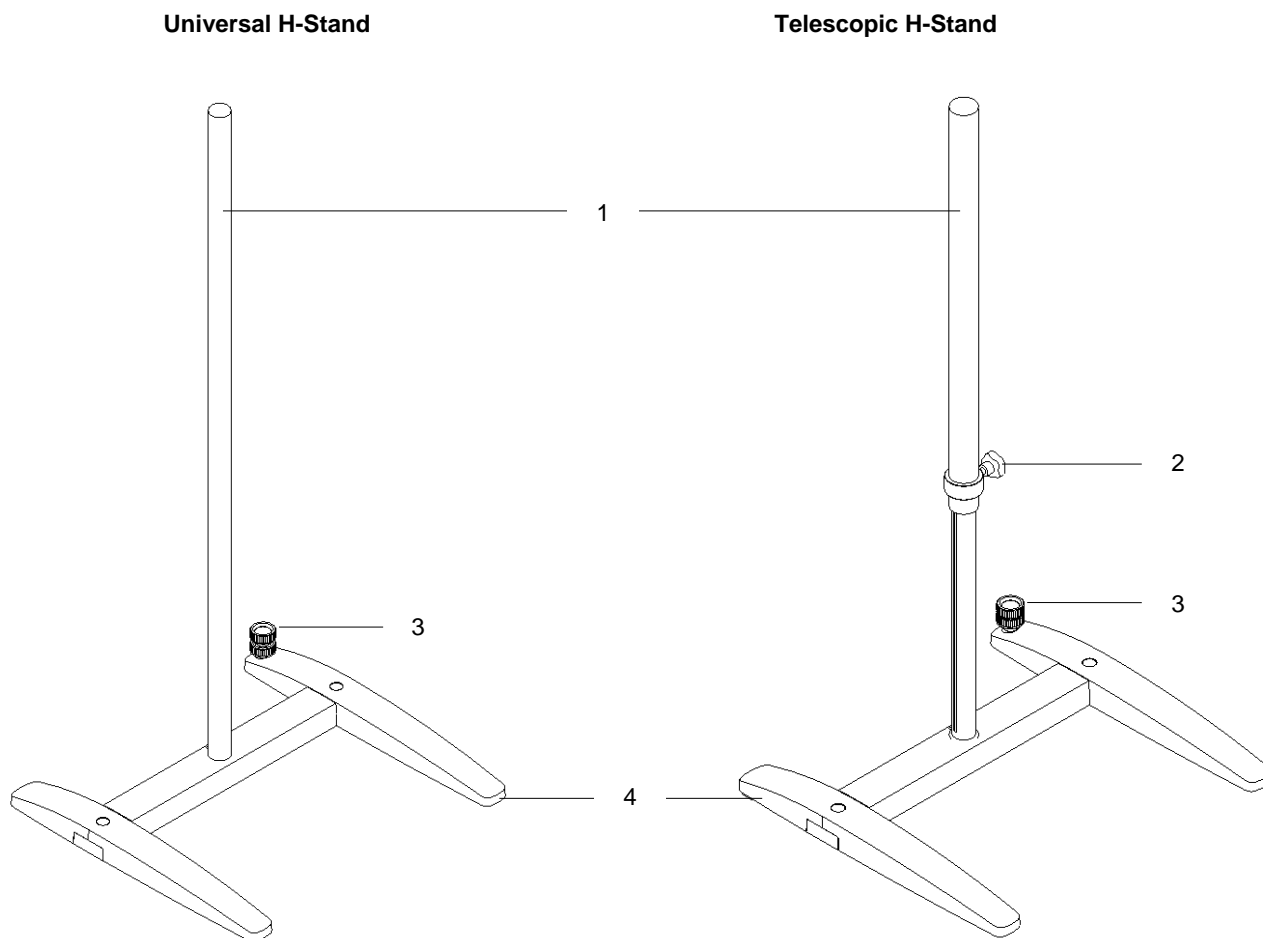
The H-Stand allow the correct fixing of overhead stirrers that require absolute stability during operation by properly using the double clamp.

Telescopic H-Stand allows the lifting of equipment in use with extreme ease and without effort by the user.

Fig. 1

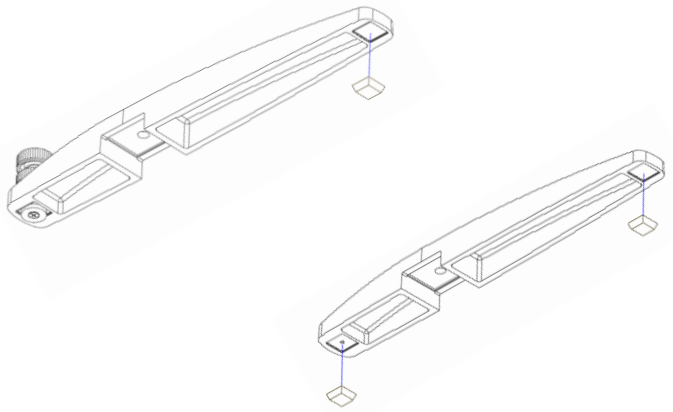
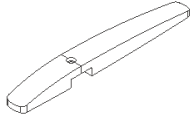
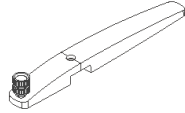
- 1 Universal Shaft  
Telescopic Shaft
- 2 Knob with screw
- 3 Adjusting screw with counter nut
- 4 Stand base

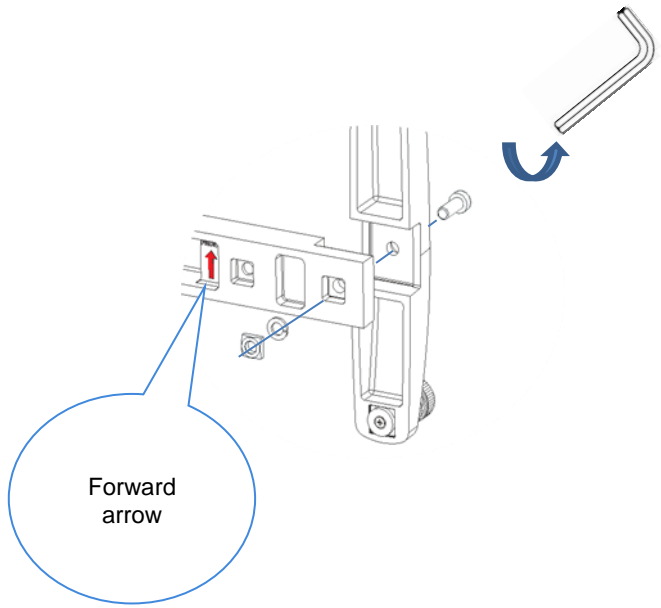
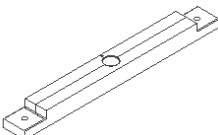
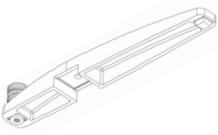
## 2. Assembly

- Unpacking
  - Check the integrity of device after the unpacking
- The box include
  - Universal H-Stand or Telescopic H-Stand
  - Instruction manual
  - Hexagonal key

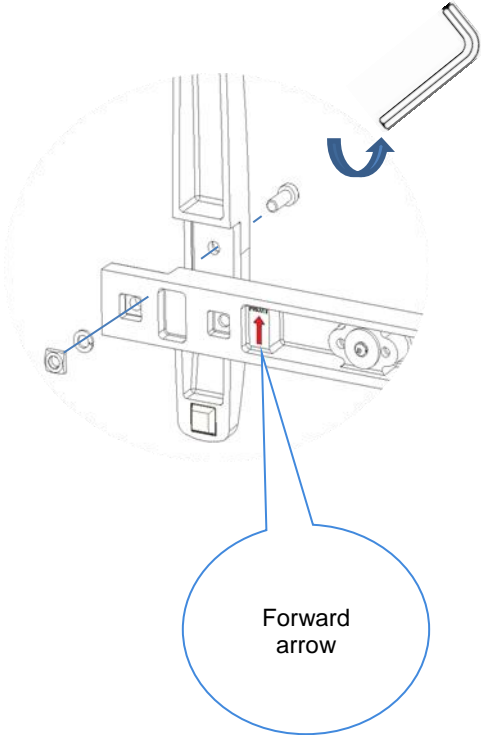
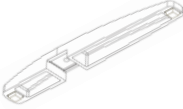
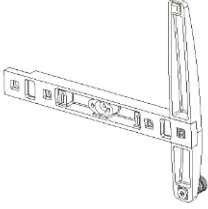
**STEP 1**

n° pieces	Description	Image part	Assembly
3	Adhesive feet		
1	Lateral H-Stand		
1	Lateral H-Stand with screw		

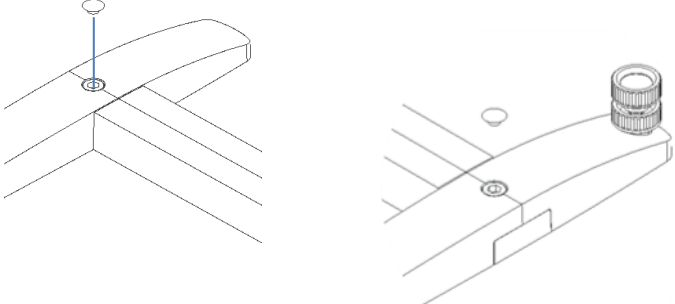
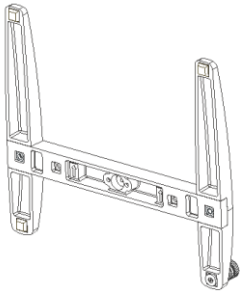
**STEP 2**

n° pieces	Description	Image part	Assembly
1	Screw M10x25 mm		
1	Nut M10		
1	Washer M10		
1	Basis H-Stand		
1	Lateral H-Stand With screw		
1	Hexagonal key 8mm		

**STEP 3**

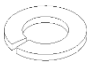
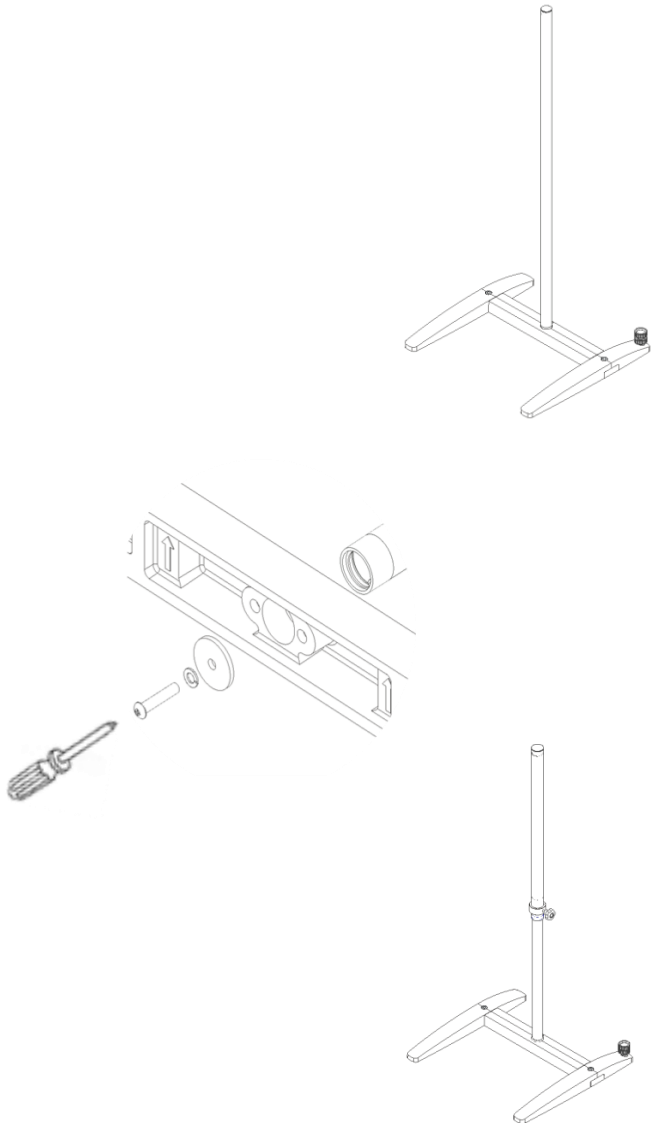
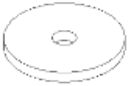
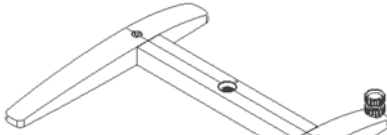
n° pieces	Description	Image part	Assembly
1	Screw M10x25 mm		
1	Nut M10		
1	Washer M10		
1	Lateral H-Stand with feet		
1	Pre- assembled H-Stand (Step 2)		
1	Hexagonal key 8mm		

**STEP 4**

n° pieces	Description	Image part	Assembly
2	Screw cover		
1	Pre- assembled H-Stand (Step 3)		

**STEP 5**

n° pieces	Description	Image part	Assembly
1	Washer M6		
1	Screw M6x30		
1	Washer M10		
1	Arm for Universal H-Stand		
1	Arm for Tescopic H-Stand		
1	Pre- assembled H-Stand (Step 4)		
1	Screw Driver NOT SUPPLIED		

**3. Installation**

- Follow this safety instruction described in this manual.
- Install the device on a flat, stable, clean, non slip, dry and fire-resistant surface.
  - With the adjustment knob with screw (Pos. 3 fig. 1) it is possible to level on an uneven surface.
  - With the lock nut it is possible to lock the adjustment knob with screw.
- Prior to each use, always check the device is in good operating condition.
- Use Double clamp properly to fix the stirrer correctly.
- Use Double clamp and Ribbon clamp properly to fix the container (Beaker with sample) correctly.
- Check the correct position of stirrer to avoid the danger of tipping over.
- Respect the max permissible load as described in the "Technical data" in this manual.

## 4. Maintenance

The Universal H-Stand and Telescopic H-Stand are maintenance free.

### Cleaning

- Cleaning should be performed with a soft cloth dampened with water containing mild surfactant. Any detergents must be approved by OHAUS.
- Wear protective gloves during cleaning of device

### Repair

If you need to contact an Authorized OHAUS Service Agent, please visit our website [www.ohaus.com](http://www.ohaus.com) to locate the Ohaus office nearest you. An Ohaus Product Service Specialist will be available to assist you.

## 5. Technical data

		Universal H-Stand 30586771	Telescopic H-Stand 30586772
Stand rod	mm	-	400
Max load	Kg		10
Arm (diameter)	mm	28	34
Dimensions (w x d)	mm	462 X 430	
Height	mm	1015	615 - 1015
Weight	Kg	4,1	4,6

## 6. Warranty

OHAUS products are warranted against defects in materials and workmanship from the date of delivery through the duration of the warranty period. During the warranty period OHAUS will repair, or, at its option, replace any component(s) that proves to be defective at no charge, provided that the product is returned, freight prepaid, to OHAUS.

This warranty does not apply if the product has been damaged by accident or misuse, exposed to radioactive or corrosive materials, has foreign material penetrating to the inside of the product, or as a result of service or modification by other than OHAUS. In lieu of a properly returned warranty registration card, the warranty period shall begin on the date of shipment to the authorized dealer. No other express or implied warranty is given by OHAUS Corporation. OHAUS Corporation shall not be liable for any consequential damages.

As warranty legislation differs from state to state and country to country, please contact OHAUS or your local OHAUS dealer for further details.

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P/N 30586845A ©2020 Ohaus Corporation, all rights reserved / todos los derechos reservados / tous droits réservés / alle Rechte vorbehalten / tutti i diritti riservati
Nous, OHAUS Corporation, déclarons sous notre responsabilité que le produit est conforme aux normes suivantes:
Ohaus Corporation
Ohaus Europe GmbH